


Instructions for Use: Pocket Microscope

Brand Name of Product	Pocket Microscope 20x–40x
Generic Name of Product	Pocket Microscope
Product Code Number(s)	MG10081-8
Purpose of Product	Ideal for onsite detailed examination of small areas and objects. Useful for quality control inspection.
Range of Applications for Product	For use with any products in need of visual inspection after cleaning.
Key Specifications of Product	<ul style="list-style-type: none"> Built in light source. Microscope stand allows ambient light to the object under examination. Camera battery included (Three [3] AG12 batteries).

Shipping & Storage	
Shipping Conditions & Requirements	N/A
Storage Conditions	N/A
Packaging Conditions	N/A
Shelf Life	N/A

Instructions for Using Product	
Description of Use(s)	Use during visual inspection of products.
Preparation for Use	Remove the Battery Insulate by pulling it outward.
Diagrams (drawings, pictures)	
Steps for Use of Product	<ul style="list-style-type: none"> Observing with Specimen Stage <ol style="list-style-type: none"> Place the object of observation onto the Specimen Slide and place a Specimen Slide Cover on top of the object. Insert the Specimen Slide into the Specimen Stage. Place the Microscope into the Specimen Stage and press the LED On/Off Button and observe through the Eyepiece. Slide the zoom lever to adjust the desired degree of magnification (from 20x to 40x). Turn the focus wheel gradually until the image is clear and sharp. Observing without Specimen Stage <ol style="list-style-type: none"> Place the Pocket Microscope directly on the top of the object of observation and press the LED On/Off Button. Slide the zoom lever to adjust the desired degree of magnification (from 20x to 40x). Turn the focus wheel gradually until the image is clear and sharp.
Interpretation of Results	N/A
Contraindications of Test Results	N/A
Documentation	N/A
Special Warnings and Cautions	<ul style="list-style-type: none"> If the LED light becomes dim, it means it is time to replace the batteries. <ol style="list-style-type: none"> Unscrew the battery door screw with a screwdriver and remove the battery door in the direction shown by the arrow. Remove the old battery and insert three (3) pcs of fresh AG12 button cells into the battery compartment in the correct polarity.

	3. Close the battery door and tighten the screw with a screwdriver.
Disposal	N/A

Reprocessing Instructions	
Point of Use	N/A
Preparation for Decontamination	N/A
Disassembly Instruction	N/A
Cleaning – Manual	N/A
Cleaning – Automated	N/A
Disinfection	N/A
Drying	N/A
Maintenance, Inspection, and Testing	N/A
Reassembly Instructions	N/A
Packaging	N/A
Sterilization	N/A
Storage	N/A
Additional Information	N/A

Related Healthmark Products	N/A
Other Product Support Documents	N/A
Reference Documents	ProSys™ Brochure, ProSys™ Price List
Customer Service Contact	Healthmark Industries Company, Inc. 18600 Malyn Blvd. Fraser, MI 48026 1-586-774-7600 healthmark@hmark.com hmark.com